

Critical-cleaning detergents for laboratory, healthcare and industrial applications

30 Glenn Street
White Plains NY 10603 USA
Tel. 914.948.4040
Fax. 914.948.4088

**24 Hour Emergency Number
(CHEM-TEL) (800) 255-3924 in
U.S.**

(e-mail) cleaning@alconox.com

(URL) <http://www.alconox.com>

- distributors
- technical information
- free samples
- new developments



Citrajel®

Low-foaming Liquid Acid Cleaner

- Ideal for use as acid wash or rinse
- Concentrated to save you money
- Phosphate free, biodegradable and readily disposable
- Low foaming acid cleaner and rinse for high performance results from your cleaning machinery
- Use to pass your cleaning validation tests for lab accreditation and plant inspection approval
- Combines acid and detergent ingredients to remove metal oxides, salts and inorganic residues

Used to clean: Laboratory ware, sampling apparatus, dairy equipment optical parts, electronic components, pharmaceutical apparatus, cosmetics manufacturing equipment, metal forgings, stampings and castings, industrial parts, pipes, tanks, boilers and reactors. USDA compliant for use in federally inspected meat and poultry plants. Passes inhibitory residue test for water analysis. FDA certified.

Used to remove: Scale, metal oxides, metal complexes, trace inorganics, milkstone, soil, grit, buffing compound, grime, grease, fats, oils, particulates, amphoteric peptides, deposits, chemical and solvents.

Surfaces cleaned: Corrosion inhibited formulation recommended for glass, metal, stainless steel, porcelain, ceramic, plastic, rubber and fiberglass. Can be used on soft metals such as copper, bronze, aluminum, zinc and magnesium if rinsed promptly. Corrosion testing may be advisable.

Cleaning method: : Laboratory warewashers, dishwashers, washer-sterilizers, washer-sanitizers, cart washers, spray washers, electrocleaning, ultrasonic cleaners flow through clean-in-place and manual cleaning if proper skin and eye protection is used.

Directions: Make a fresh 1 - 2% solution (3 - 5 Tbsp. per gal., 1 1/4 - 2 1/2 oz. per gal. or 10 - 20 ml per liter) in cold, warm or hot water. If available, use hot water. For difficult soils, raise water temperature and use more detergent. Clean by soak, circulate, wipe or ultrasonic method. Wear protective gloves and eyewear. RINSE THOROUGHLY--preferably with running water. For critical cleaning, do final or all rinsing in distilled, deionized or purified water. For food contact surfaces, rinse with potable water. Used on a wide range of glass, ceramic, plastic and metal surfaces. Corrosion testing may be advisable.

Available in Convenient Sizes

	Alconox Cat. #
Case 4 x 1 Gallons	2001
15 Gal. Drum	2015
30 Gal. Drum	2030
55 Gal. Drum	2055
1 gal. makes 50 gal. cleaning solution	



Citrajel is available from leading laboratory, hospital, clinical and industrial suppliers. To find a distributor for Alconox, Inc. detergents, visit "Find Dealer" at the website. To request FREE samples, visit Sample Request at www.alconox.com, write or call Alconox, Inc. today.

- distributors
- technical information
- free samples
- new developments

PHYSICAL DATA

TYPICAL VALUE

pH of 1% solution	2.5
Specific gravity (g/ml)	1.10
Density (lbs./gal.)	9.15
Flash Point (degrees F):	None
Phosphate Content (as Phosphorus)	0%
Other Phosphorous compounds (as Phosphorous)	less than 1%
Organic Carbon (% calculated w/w)	14%
Fragrance Content	0%
Surface Tension 1% Solution (Dyne/cm)	32
Color: Pale Yellow	
Form: Liquid	
Solubility in Water: Completely soluble in all proportions	
Hard Water Effectiveness: Highly Effective	
Biodegradability: Biodegradable	
Foam Tendency: Low to Non- Foaming	
Shelf Life: Two years from the date of manufacture	

Chemical Description

Citrajel contains a blend of organic acids and surfactants. Citrajel is phosphate free.

Cleaning Validation Methods:

Test a parameter of rinse water before and after rinsing the cleaned surface, or test the clean surface. No significant change in the parameter indicates no detectable residue. Parameters measured include: pH, conductivity, total organic carbon, surface tension, and surface analysis.

To get a bulletin on cleaning validation method references for Alconox, Inc. detergents, go to the Pharmaceutical Cleaning Validation References page in the TECH INFO section of our website at www.alconox.com.

Health Safety Information:

OSHA Hazardous Ingredients: None
RCRA Hazard Class: Non-hazardous
EPA Priority Pollutants: None
DOT Hazard Class: Non-hazardous
Flammability: Non-flammable
Oral Toxicity: LD₅₀ > 5000 mg/kg oral rat. No ingredient defined as an oral toxicant by OSHA
Eye Irritation: Serious potential for eye irritation if not rinsed
Inhalation Toxicity: Non-irritating
Latex Content: no latex ingredients or latex in packaging
Carcinogenicity:
NTP = No IARC = No OSHA = No
VOC Content: 0%

All ingredients in Citrajel are listed in TSCA inventory.

Precautions:

No special precautions other than good industrial hygiene and safety practices employed with any industrial chemical (see Directions). A Material Safety Data Sheet is available at www.alconox.com or by calling fax-on-demand at 914-948-4040 and following the prompts. Use fax document #0320.

Contact Alconox, Inc. for purchase specifications. Typical data is not a specification.

While the information in this report should not be considered to be a product warranty, we urge you to investigate, test and verify the suitability of Alconox detergents for your specific application. We, of course, can not give permission to use, or recommend the use of, our detergents where they infringe patents. No representation or warranty is made as to the safety of products or materials mentioned under the Federal Food Additives Amendment of 1958.